

Injectable Grade of Sucrose

What is sucrose ?


- A disaccharide consisting of glucose and fructose joined by α -1,2 glycosidic bonds, it is mainly refined from sugar beet and sugarcane.
- Glucose's aldehyde group and fructose's ketone group form a glycosidic bond, so it does not have reducing properties and is considered a non-reducing sugar.
- A colorless crystal.
- Extremely soluble in water.
- Slightly soluble in methanol and insoluble in ethanol (95%).



SUCROSE SG

Sucrose structure

Chemical Formula : $C_{12}H_{22}O_{11}$
MW : 342.30
CAS RN® : 57-50-1


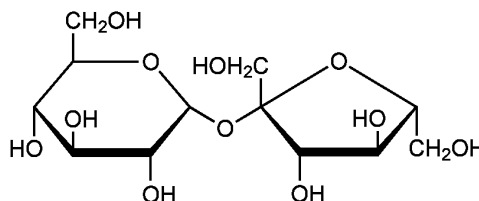


Applications and references

- Improved stability of mRNA LNP formulations against freeze-drying.
 - Lamoot A., *et al.* Biomater Sci. 11(12), p4327-4334 (2023)
Ex.: Lipid Nano Particle (LNP) formulation
 - Importance of selecting pharmaceutical additive sucrose with low nanoparticle impurities (NPI).
 - Weinbuch D., *et al.* Pharm. Res. 32, p2419-2427 (2015)
 - Weinbuch D., *et al.* Pharm. Res. 34, p2910-2921 (2017)
Ex.: Therapeutic antibody
 - Used as a freezing/cryoprotectant for liposomes.
 - Guimarães D., *et al.* Front. Bioeng. Biotechnol. 7, Article 424 (2019)
Ex.: Freezing/cryoprotection of liposomes containing tamoxifen.
- ◆ IUPAC name : β -D-Fructofuranosyl α -D-glucopyranoside

The information provided herein is intended only for reference purposes. It is the customer's responsibility to determine that the ingredients meet all legal requirements in the country where they are used, and that they do not infringe on any third party patents. Unauthorized reproduction of this brochure is prohibited.

Application study for lipid nano particles (LNP)

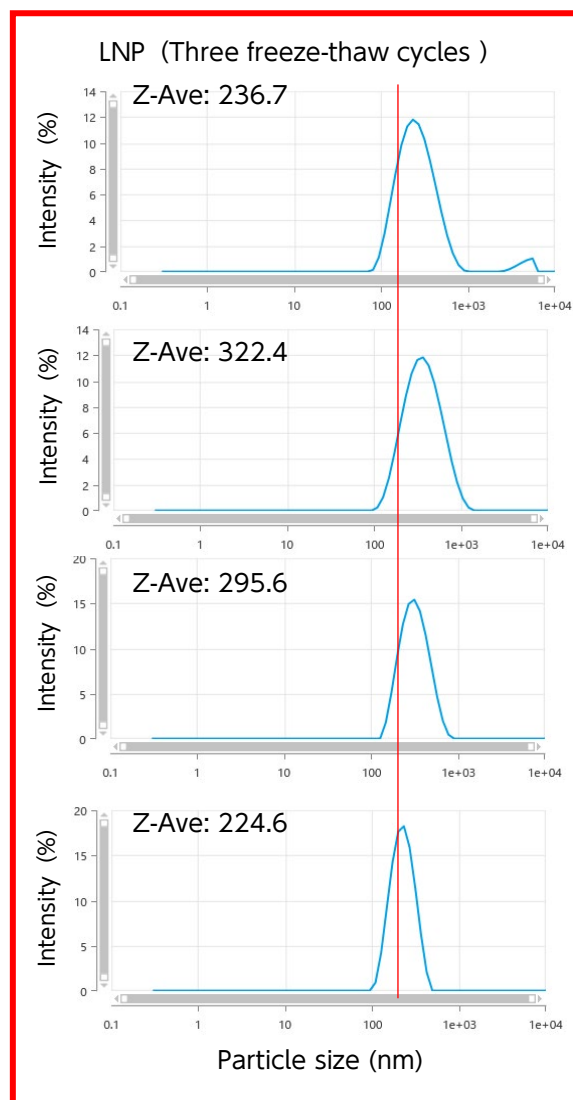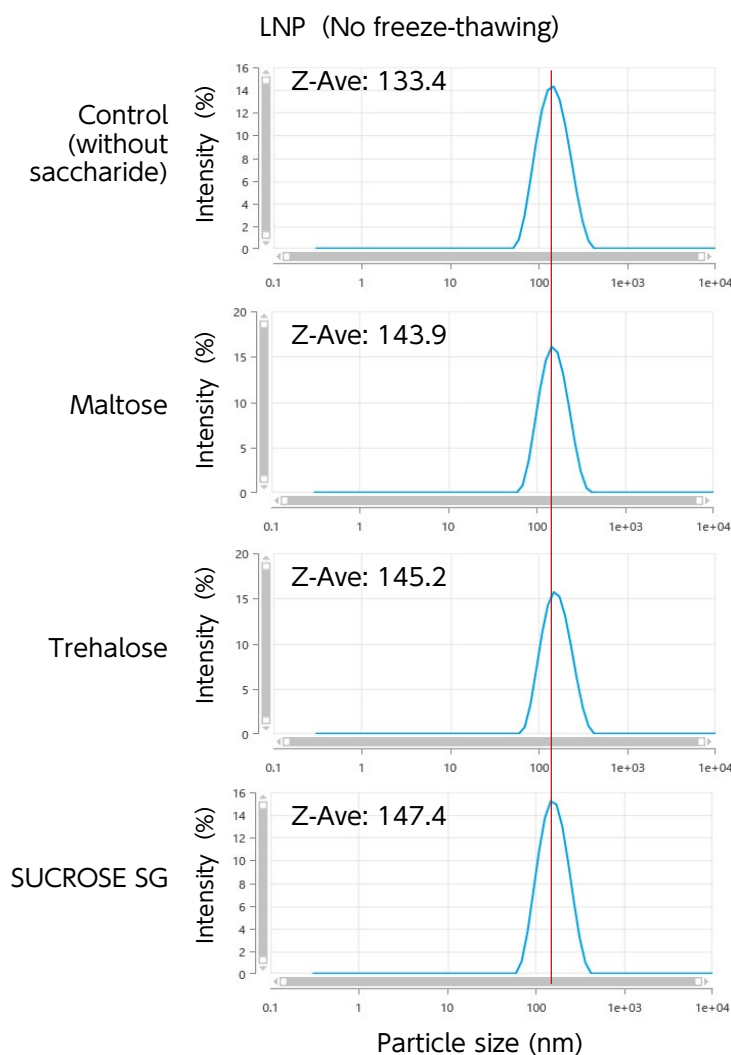

【Materials and Methods】

Prepared using Cayman's LNP preparation kit; LNP compositions were prepared by replacing sucrose in the Comirnaty® composition with the respective carbohydrate and subjected to freeze-thaw testing.

【Results】

After the freeze-thawing (-20°C, 1 hr ⇔ R.T., 0.5 hr), the addition of SUCROSE SG resulted in the least change in particle size. A trend towards larger particle size of LNP was observed under the others.

*LNP stabilization by freeze-thaw cycles.

(SUCROSE SG > Trehalose > Maltose > without saccharide)

MANUFACTURER : **Nagase Viita Co., Ltd.**
CONTACT : **Nagase & Co., Ltd.**

Life & Healthcare Products Dept. Pharma-Medical Div.
TEL: +81-3 (3665) 3333 (TOKYO JAPAN) TEL: +81-6 (6535) 2327 (OSAKA JAPAN)
E-mail: dnfct@ex.nagase.co.jp

⚠ The information provided herein is intended only for reference purposes. It is the customer's responsibility to determine that the ingredients meet all legal requirements in the country where they are used, and that they do not infringe on any third party patents. Unauthorized reproduction of this brochure is prohibited.